

GENEL[®]

Gas Generator Technologies

DESIGN AND
PRODUCTION OF
LABORATORY
GAS
GENERATORS

N₂ H₂ O₂ ZERO AIR



**ZERO AIR
GENERATORS**

ZA series

ZA Series Zero Air

The CINEL ZA series Air Generators produce laboratory grade purified and hydrocarbon free air from an existing in house oil free compressed air supply, eliminating the need for inconvenient high-pressure gas cylinders. The CINEL ZA Series Zero Air Generators remove HC pollutants to less than 0.1 ppm, CO to less than 0.1 ppm and all kinds of particles. The system is designed to be a complete turnkey, is easy to install, and requires only minimal annual maintenance. The pretreated air passes through a coalescing filter to eliminate oil, water and the particles higher than 5 microns. The hydrocarbons including methane are removed by oxidation on a platinum/palladium catalyst heated at high temperature. A particle filter at the exit of the catalyst, makes it possible to eliminate the residual particles from outgoing flow.

Application

GC-FID
GC-FPD, GC-NPD



Technical data

	ZA 1500	ZA 3000	ZA 6000
FLOW (l/min)	1.5	3	6
HC & CO out	< 0.1 ppm	< 0.1 ppm	< 0.1 ppm
MAX CO in	50 ppm	50 ppm	50 ppm
MAX HC in	100 ppm	100 ppm	100 ppm
MAX TEMP. in	40 °C	40 °C	40 °C
PRESSURE in	4.5 - 10 bar	4.5 - 10 bar	4.5 - 10 bar
PRESSURE drop	< 1 bar	< 1 bar	< 1 bar
STABILITY (min)	40	40	40
In / out CONNECTIONS	1/4 - 1/8 (NPT)	1/4 - 1/8 (NPT)	1/4 - 1/8 (NPT)
POWER	220 V / 50 Hz or 110 V / 60 Hz	220 V / 50 Hz or 110 V / 60 Hz	220 V / 50 Hz or 110 V / 60 Hz
WEIGHT (kg)	18	18	18
DIMENSIONS (cm)	28x43x17H	28x43x17H	28x43x17H
WORKING TEMP.	Amb. +20°C	Amb. +20°C	Amb. +20°C

OTHER PRODUCTS

NITROGEN GENERATORS **N₂**

HYDROGEN GENERATORS **H₂**

OXYGEN GENERATORS **O₂**

CINEL Strumenti Scientifici s.r.l.
via dell'Artigianato, 14-14/A, 35010 Vigonza (Padova) - Italy
tel. +39 049 725022 fax +39 049 8931881 e-Mail info@cinel-gas.com
www.cinel.com - www.cinel-gas.com



SINCERT

Azienda con sistema di qualità certificato ISO 9001:2008

Wyłączny dystrybutor: Genore chromatografia, e-mail: info@genore.pl, www.genore.pl, tel. 22 40 107 342